

EasyDirect Pipette Asset Management

Getting Started is Easy

If you're already tracking your pipette inventory with a spreadsheet, your current .xls, .csv and other formatted files upload directly into EasyDirect. The software is designed to optimize and adapt to the way you're currently managing your pipettes.



Version	Starter*	Site	Unlimited
Catalog number	30475911	30475913	30475915
Pipettes	200	1500	Unlimited
Connected SmartStands	5	200	Unlimited
Calibration certificate upload	Manual Upload	Automatic**	Automatic**
Installation type	Single client	Single client/server	Single client/server

*Included with SmartStand **Check for local availability

	Computer	Minimum	Optimum
Database Server	CPU	2 GHz Dual Core	3 GHz Quad Core
	RAM	4 GB	16 GB
	Display Resolution	1280 x 800	1280 x 1024
	Hard Disk	30 GB free space	100 GB free space
	Operating System	Windows 7 SP1	Windows Server 2016
	SQL Server	MS SQL Server 2008 R2 SP3	MS SQL Server 2016
EasyDirect PAM Business Server	CPU	2 GHz Dual Core	3 GHz Quad Core
	RAM	4 GB	8 GB
	Display Resolution	1280 x 800	1280 x 1024
	Hard Disk	30 GB free space	10 GB free space
	Operating System	Windows 7 SP1	Windows Server 2016
	SQL Server	MS SQL Server 2008 R2 SP3	MS SQL Server 2016
EasyDirect Client	CPU	2 GHz Dual Core	3 GHz Quad Core
	RAM	4 GB	8 GB
	Display Resolution	1280 x 800	1920 x 1080



See it in action!
mf.com/EDPAM-demo

Asset Management



Pipette Asset Management

Comprehensive Tracking

Optimize Calibration Scheduling

Scalable – Lab to Enterprise

Be Audit-Ready... Always!

www.mt.com/easydirectpam

For more information

Save Time, Money and Frustration

Pipette Management Simplified

METTLER TOLEDO Group
Laboratory Division
Local contact: www.mt.com/contacts

Subject to technical changes
© 08/2018 METTLER TOLEDO. All rights reserved.

17700929_EN Rev A

Distributed By:



Adelab Scientific
36 Holland Street
Thebarton SA 5031
Ph 08 8234 7955
Fax 08 8234 7897
info@adelab.com.au
www.adelab.com.au


METTLER TOLEDO

With EasyDirect, You're in Control

Save Time and Simplify Pipette Management

With EasyDirect Pipette Asset Management (EDPAM) and SmartStand, you have a powerful pipette tracking and calibration management system. Using the RFID chip embedded in every Rainin XLS-brand pipette, EasyDirect tracks every pipette's usage, location and service data automatically.

- Manage an unlimited number of pipettes
- Track all pipettes automatically
- Simplify calibration management and repairs

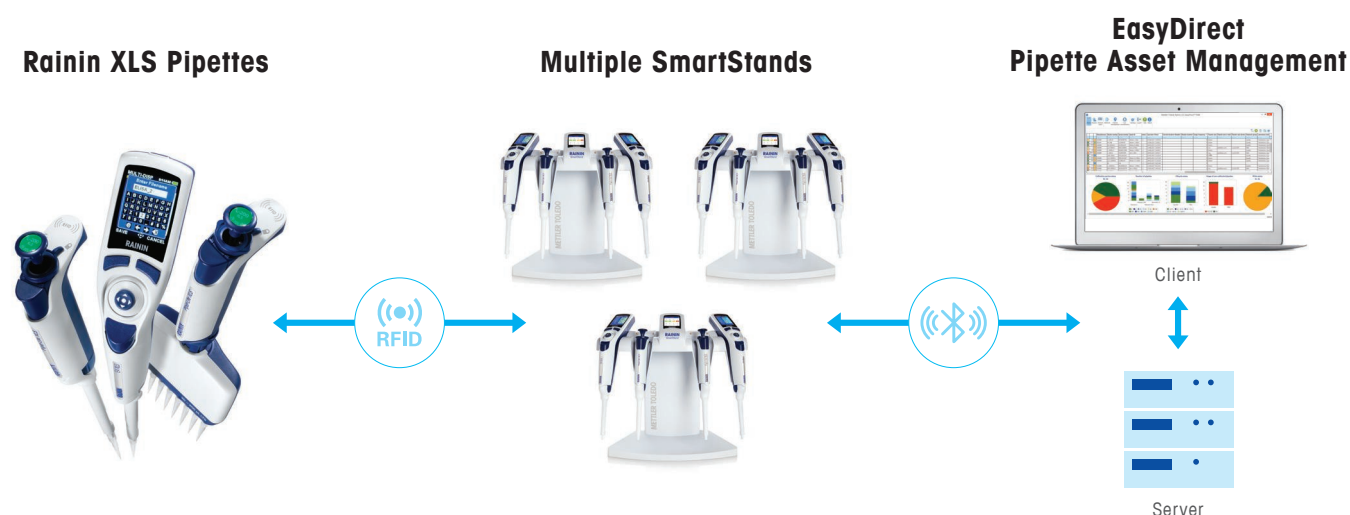


A simple, unified system for tracking and managing pipettes (all manufacturers)

Place any Rainin XLS-brand pipette on a SmartStand to instantly see the pipette's calibration status and service details right at the bench. EasyDirect connects to SmartStand via Bluetooth or USB drive to automatically transfer serial numbers, service and other data to the EasyDirect database.

EasyDirect maintains a complete history of your entire pipette inventory including calibration certificates and verification reports. This makes it easy to track the status, location, calibration and service schedule for any pipette.

That's pipette management...simplified!

- Assign locations, users and groups to any pipette
- See all pipette details on a single screen
- Search and filter – create custom reports and schedules
- See the current calibration/maintenance status of all, some or even just one pipette
- Minimize disruptions – EasyDirect helps plan and manage calibration
- Setup automatic email service reminders for end-users or support groups

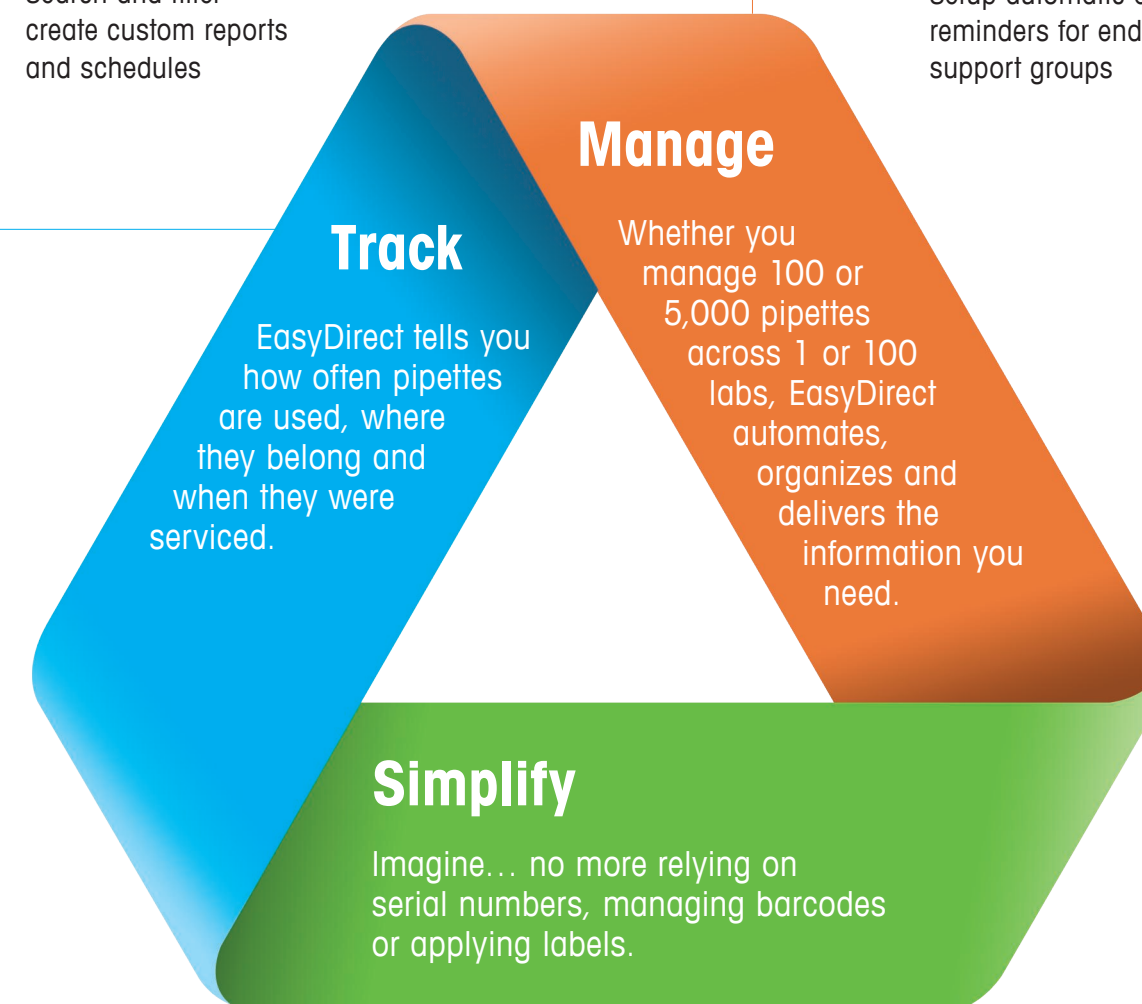

- EasyDirect uses SmartStand to automatically transfer serial numbers, service data and other information to and from your pipettes
- EasyDirect maintains a comprehensive history for each pipette, automatically assigning service and verification records.
- Let EasyDirect update the last and next calibration dates automatically for you after each service, saving you valuable time.

Trust the Service Experts

Minimize Risk and Maximize Reproducibility

The METTLER TOLEDO difference: controlled laboratory environments and manufacturer authorized spare parts, certified pipette technicians and specialized calibration software.



Uptime
Our global network of factory-trained technicians and ISO/IEC 17025 certified service labs enables us to respond quickly to all repair and calibration needs.



Performance
Environmentally controlled labs and original manufacturer spare parts ensure that instruments are always returned to the highest accuracy and precision.



Compliance
With specialized calibration software and ISO/IEC 17025 accredited labs, METTLER TOLEDO is a partner you can trust to help maintain ISO 8655 and FDA 21 CFR part 11 standards.



Expertise
Our certified pipette service technicians perform more than 10 million readings each year, giving us extraordinary data analytics for understanding how all brands of pipettes perform over time.

METTLER TOLEDO operates the world's most extensive network of ISO/IEC 17025 accredited pipette calibration laboratories.

- Comprehensive offering**
- Calibration
 - Preventive maintenance
 - Service repair
 - Calibration certificates

Global network
With a global network of ISO/IEC 17025 accredited laboratories and one of the largest factory-trained support and service forces, we can provide you with application support and technical assistance anytime, anywhere.

► Request Rainin Service today! www.mt.com/RaininService

Routine Pipette Checks Simplified

Periodic Checks Assure Performance

Periodically assessing a pipette is performing within tolerances ensures data integrity and reduces risk.

You can turn a conventional balance into a pipette check station by simply mounting an evaporation trap onto the weighing pan. The onboard application on the METTLER TOLEDO XPE and XSE balance guides users through the testing procedure and records each pipette’s test results, which it sends electronically to the EasyDirect PAM software.



How EDPAM simplifies your pipette management

